

## TM 232 Bearing friction



### Learning objectives/experiments

- determine the frictional moment in slide bearings with various friction pairs
- determine the frictional moment of a rolling bearing
- comparison of slide and rolling bearings
- basic experiments on rotational dynamics

#### Description

- friction in slide bearings and roller bearings
- interchangeable bearing shells for slide bearings of different materials

Bearings can be classified as slide bearings or roller bearings depending on the nature of their motion. In slide bearings, there is a sliding motion between the bearing and the supported component. In roller bearings, there is both sliding motion and rolling motion between rolling bodies and the supported component. Sliding motion is undesirable in roller bearings, e.g. between rolling body and cage. The bearing force is transferred in the roller bearing by means of rolling. In both types of bearings—the slide bearing and the roller bearing—frictional forces occur during operation, which oppose resistance to the movement.


The TM 232 experimental unit allows investigations of friction on slide bearings with different bearing shells and on roller bearings. A shaft with a cable drum and flywheel is mounted on a base plate. The weight of the heavy flywheel generates bearing forces. A moment is applied by means of weights, which is equal to the friction moment at the start of the rotation. Replaceable bearing shells are used as slide bearings. The coefficients of friction are determined in experiments. Bearing shells made of different materials are included in the scope of delivery in order to study different friction pairings. The bearing friction is very low when using the roller bearing. In this case, the flywheel can be used for basic experiments on rotational dynamics.

The experimental unit is designed to be fixed to a wall. The parts of the experiment are clearly laid out and securely housed in a storage system.


G.U.N.T. Gerätebau GmbH, Hanskampring 15-17, D-22885 Barsbüttel, Telefon (040) 67 08 54-0, Fax (040) 67 08 54-42, Email sales@gunt.de, Web www.gunt.de We reserve the right to modify our products without any notifications. Page 1/2 - 03.2017




# **TM 232** Bearing friction

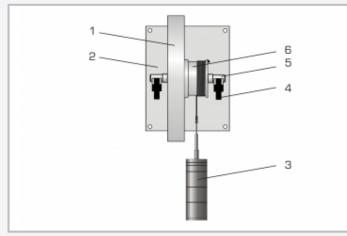

1 flywheel, 2 base plate, 3 weight, 4 bearing seat, 5 shaft, 6 cable drum

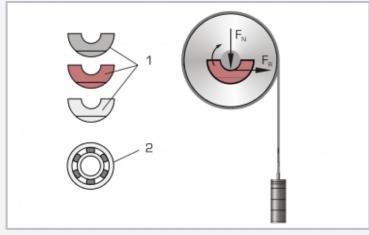

1 interchangeable bearing shells of cast iron, red bronze and plastic (PTFE), 2 roller bearing;  $F_N$  normal force,  $F_R$  frictional force

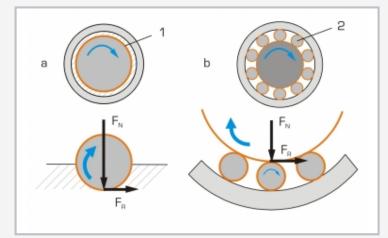

a dynamic friction in the slide bearing, 1 sliding surface

b dynamic friction and rolling friction in the rolling bearing, 2 roller bearing;  $F_{\rm N}$  normal force,  $F_{\rm R}$  frictional force

## Specification

- [1] comparison of dynamic friction and rolling friction
- [2] experiments on rotational dynamics are possible
- [3] bearing shells of different materials as slide bear-
- ings [4] steel flywheel, galvanised
- [5] drive via cable drum and weights
- [6] storage system for parts
- [7] bracket for wall mounting

### Technical data

Bearing shells as slide bearing, half-shells ■ GG-25

- red bronze
- PTFE (Teflon)

Deep-groove ball bearing ■ type 6203

Shaft bearing journal ■ Ø=17mm

Flywheel

- ∎ Ø=300mm
- weight: 22,2kg

Weights

- 1x 1N (hanger)
- 5x 1N
- 1x 2N
- 3x 5N

Base plate

LxW: 250x200mm

LxWxH: 200x330x300mm Weight: approx. 30kg LxWxH: 290x140x130mm (storage system)

Scope of delivery

- 1 experimental unit
- 6 bearing shells
- 2 roller bearings
- 1 set of weights
- 1 storage system
- 1 set of instructional material