

# WP 310

Materials testing, 50kN



The illustration shows WP 310 together with the accessory WP 310.05.

#### Description

- hydraulically operated trainer, based on industrial standards
- direct generation of tensile and compressive forces
- extensive accessories for experiments from destructive materials testing

A classic discipline of materials testing is the destructive testing method, in which specimens are mechanically tested to failure. The materials test provides data for hardness, rigidity and strength in a reproducible and precisely quantified manner.

The WP 310 unit, in conjunction with the accessories, offers experiments from destructive materials testing.

The clean layout and simple operation mean the experimental sequence can be observed in all details and phases. The power of the trainer allows tests to be performed on an industrial scale. Material specification data and laws can be verified using self-determined measured values.

The vertical, hydraulically operated trainer with direct force generation can produce both tensile and compressive forces. The height of the lower crossmember can be adjusted for coarse adjustment. Cylindrical receptacles on the cross-members allow for easy exchange of accessories. The extensive accessories provide tensile and compression tests, Brinell hardness tests, bending, shear and cupping tests. Plate and coil springs can also be tested. The test load and elongation of the specimen are measured by sensors and are displayed.

The measured values are transmitted directly to a PC where they can be analysed using the software included.

## Learning objectives/experiments

- together with the accessories
  - tensile test
  - compression test
  - Brinell hardness test
  - bending test
  - ▶ shear test
  - cupping test
  - spring testing


G.U.N.T. Gerätebau GmbH, Hanskampring 15-17, D-22885 Barsbüttel, Telefon (040) 67 08 54-0, Fax (040) 67 08 54-42, Email sales@gunt.de, Web www.gunt.de We reserve the right to modify our products without any notifications. Page 1/3 - 08.2017



# WP 310 Materials testing, 50kN



1 hydraulic cylinder for generating tensile and compressive forces, 2 operating area with the accessory WP 310.05, 3 force sensor, 4 adjustable height lower cross-member with lock, 5 displays and controls, 6 displacement sensor

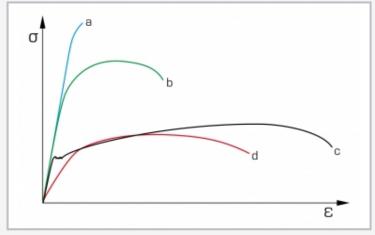

Stress-strain diagram for various materials: a hardened steel, b tempered steel, c annealed steel, d alloyed aluminium



Software screenshot: Brinell hardness test

### Specification

- hydraulically operated trainer for materials testing, based on industrial standards
- [2] generation of tensile and compressive forces
- [3] adjustable test load and travel velocity
- [4] generation of test load via gear pump and doubleacting hydraulic cylinder
- [5] force measurement via a strain-gauge full bridge with acoustic overload signal, max. overload 150%
- [6] displacement measurement via linear potentiometer
- [7] LED displays for force and displacement with tare and maximum-value storage
- [8] GUNT software for data acquisition via USB under Windows 7, 8.1, 10
- [9] wide range of accessories available

#### **Technical data**

Operating area, WxH: 300x925mm Hydraulic generation of the test load ■ test load: 0...50kN

- LEST IDAU. U...JUKIN
- max. system pressure: 175bar
- max. piston stroke: 150mm
- traverse velocity: 0...425mm/mingear pump
  - ▶ max. flow rate: 1 cm<sup>3</sup>/revolution
  - ▶ power consumption: 0,55kW

#### Measuring ranges

- force: 0...50kN
- displacement: 0...150mm

230V, 50Hz, 1 phase 230V, 60Hz, 1 phase; 120V, 60Hz, 1 phase UL/CSA optional LxWxH: 1080x830x2300mm Weight: approx. 330kg

### Required for operation

PC with Windows recommended

#### Scope of delivery

- 1 trainer
- 1 GUNT software CD + USB cable
- 1 set of instructional material



# WP 310 Materials testing, 50kN Hydraulic universal material tester, 50kN

Optional accessories

| Tensile test              |           |   |
|---------------------------|-----------|---|
| 020.31005                 | WP 310.05 | Wedge grips for tensile specimens, round and flat   |
| 020.31012                 | WP 310.12 | Set of 10 tensile specimens F10x50, St              |
| or                        |           |   |
| 020.31006                 | WP 310.06 | Threaded grips for tensile specimens, threaded end  |
| 020.31011                 | WP 310.11 | Set of 10 tensile specimens B10x50, St              |
| or                        |           |   |
| 020.31007                 | WP 310.07 | Holder for tensile specimens, dumbbell-shaped       |
| 020.31013                 | WP 310.13 | Set of 10 tensile specimens, dumbbell-shaped, St    |
| Compressive strength test |           |   |
| 020.31004                 | WP 310.04 | Compression plates for compression tests            |
| 020.31015                 | WP 310.15 | Set of 5 compression specimens, 4x plastic, 1x wood |
| Brinell hardness testing  |           |   |
| 020.31001                 | WP 310.01 | Experimental setup for Brinell hardness tests       |
| 020.30003                 | WP 300.03 | Set of 4 hardness specimens, Al, Cu, St, CuZn       |
| 020.30031                 | WP 300.31 | Set of 4 hardness specimens, Al                     |
| 020.30032                 | WP 300.32 | Set of 4 hardness specimens, Cu                     |
| 020.30033                 | WP 300.33 | Set of 4 hardness specimens, St                     |
| 020.30034                 | WP 300.34 | Set of 4 hardness specimens, CuZn                   |
| Bending test              |           |   |
| 020.31003                 | WP 310.03 | Bending test device                                 |
| 020.31084                 | WP 310.84 | Set of 25 bending cpecimens, cast iron (EN-GJL-250) |
| Cupping test              |           |   |
| 020.31010                 | WP 310.10 | Device for cupping tests                            |
| 020.30041                 | WP 300.41 | Set of 5 cupping specimens, Al                      |
| 020.30042                 | WP 300.42 | Set of 5 cupping specimens, Cu                      |
| 020.30043                 | WP 300.43 | Set of 5 cupping specimens, St                      |
| 020.30044                 | WP 300.44 | Set of 5 cupping specimens, CuZn                    |
| Shear test                |           |   |
| 020.31002                 | WP 310.02 | Device for shear tests                              |
| 020.30052                 | WP 300.52 | Set of 5 shear specimens, Cu                        |
| Spring tests              |           |   |
| 020.31008                 | WP 310.08 | Experimental setup for spring tests, helical spring |
| 020.31009                 | WP 310.09 | Experimental setup for spring tests, disk spring    |